

Installing Windows 2000/XP 4pi FPXDriver

[\[download this section as pdf\]](#)

- Open the **Drivers** folder on the CD.
- Open the **FPXDriver** folder.
- Perform the following steps to install the Win2k-compatible FPXDriver:
 - After the Spectral Engine Hardware is installed and the computer booted, the computer will (supposedly) detect that new hardware has been installed. The Wizard will be automatically launched by Windows and display the **Found New Hardware** window.
 - Click **Next** to continue. The installer should automatically find the CD-ROM and select the correct file. If it does, you will be able to click **Next** through to the **Finish** of the procedure.
 - If the procedure does not find the correct files, tell the Wizard to **search** for a suitable driver for the device and click **Next**.
 - Uncheck all search check-boxes. Turn **on** the **Specify a Location** checkbox and hit **Next**.
 - In the subsequent search window, hit **Browse** and select the **FPXDriver** folder in the **Driver** folder on the CD-ROM.
 - The Wizard should find the FPXDrv.inf script file. Click **Open** to insert the path into the **Found New Hardware** Wizard and **OK** to confirm the selection. The **Other PCI Bridge Device** should be listed as "found." In this **Results** window press **Next** to proceed with the installation.
 - The completion window should appear. Press **Finish** to close the Wizard. No restart is necessary.
 - Install the hasp driver by running the **Hasp32.exe** installer in the **Hasp** folder, located in the Drivers folder on the CD-ROM. Don't forget to connect the hasp. Refer to the Hardware Install procedures for more information.
 - Install the FingerLakes USB driver by running the **FLIUpdate** installer in the **CCD** folder, located in the Drivers folder on the CD-ROM. Don't forget to connect the CCD camera. Refer to the Hardware Install procedures for more information.