

Installing/Updating Windows NT 4pi FPXDriver

[\[download this section as pdf\]](#)

- **Log on** to the computer as Administrator.
- Open the **Drivers** folder on the 4pi CD-ROM or software download.
- Open the **Windows NT4** folder on the 4pi CD-ROM or software download.
- Open the **FPXDriver** folder. It should look similar to that shown below:



- Run Install.BAT
- Shutdown the computer and connect the hasp if it has not been connected.
- Reboot the computer.

*Note: this procedure applies to first time installs or reinstalls. If you **update** the driver using this procedure, Revolution must not be running when the update is performed. It is of course possible to use the Device Manager to stop the driver and restart it after the update without a reboot, but this procedure is not supported.*