


The Multi-Toolbar

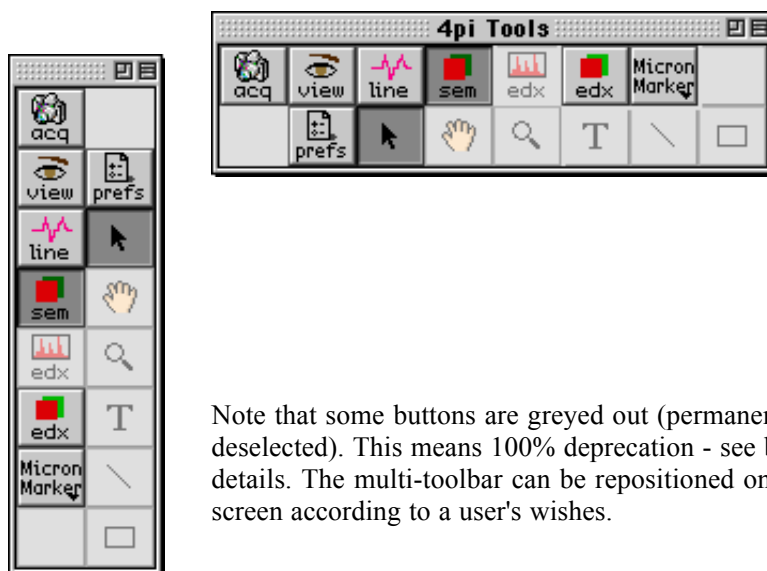



The Multi-Toolbar is a remnant of earlier versions of Revolution. As Revolution has continued development, the functionality of the toolbar has been moved to other parts of the program. This page summarizes its remaining use. The multi-toolbar will disappear completely sometime in the 1.6.x development cycle.











Select the **multi-toolbar** from the bottom of the **Mode** menu. A floating toolbar is displayed on screen as shown below. The default orientation of the toolbar is vertical. The orientation can be set with a menu in the [General Prefs](#).



Note that some buttons are greyed out (permanently deselected). This means 100% deprecation - see below for details. The multi-toolbar can be repositioned on the screen according to a user's wishes.

The function of each multi-toolbar button is described below, with links (where appropriate) to a corresponding section of the manual. *Deprecated functions may still work but are not supported.*

Button	Function	Status
	Initiate acquisition of images, x-ray maps, or spectra	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Required for x-ray mapping .
	Open preview windows of images, x-ray maps, or spectra	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Required for x-ray mapping .
	Turn linescanning on/off	Required for normal line-scanning . Required for X-ray line-scanning .
	Select SEM images function.	Deprecated. Use SEM Image Mode .

	Select EDS Spectrum for acquisition.	Deprecated. Use EDX Spectrum Mode .
	Include x-ray image maps in acquisition.	Required for x-ray mapping .
	Turn on/off the micron marker.	Deprecated for SEM Imaging (use SEM Image Mode and Micron Marker Prefs). Still used for x-ray mapping .
	Open the preference panels	Deprecated. Use Edit → Preferences... menu.
	Switch to pointer-cursor, for selecting items	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Still used for x-ray mapping .
	Switch to hand-cursor, for dragging items.	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Still used for x-ray mapping .
	Switch to magnify-cursor, for zooming items.	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Still used for x-ray mapping .
	Switch to text-cursor, for editing.	Deprecated for SEM Imaging (use SEM Image Mode). Deprecated for EDX Spectra (use EDX Spectrum Mode). Still used for x-ray mapping .
	Select line drawing cursor.	Never implemented.
	Select rectangle drawing cursor.	Never implemented.