

Removing WinRT from Windows NT

[\[download this section as pdf\]](#)

Removing WinRT from Windows NT

- **Log on** to the computer as Administrator.
- **Disable WinRT.**
 - Open the **Control Panels** folder and double-click **Devices**.
 - Select **WinRT** from the list (abort this step if WinRT is not listed).
 - Click **Startup** and click the **Disabled** radio button.
 - Click **OK** and close the **Devices** window.
 - Reboot the computer.
- **Remove WinRT.** To do this you must know the location of the **systemroot** folder. This is on the boot drive and likely is called "C:\Winnt," but may be called something else depending on your installation.
 - Open **systemroot\system32\drivers**
 - Delete the file named **WinRT.sys** (abort this step if WinRT.sys is not present)
 - Use the Run command in the START menu to run **regedit**
 - Navigate the registry tree to
HKEY_LOCAL_MACHINE→System→CurrentControlSet→Services→WinRT
 - Delete the **WinRT** folder.
 - Close **regedit**.
- Proceed to the [4pi FPXDriver installation procedure](#) appropriate for your operating system.

Removing WinRT from Windows 98SE

- **Disable WinRT.**
 - Open the **Control Panels** folder and double-click **System**.
 - Click the **Device Manager** tab and view **by type**.
 - Expand the WinRT device category (abort this step if WinRT is not listed).
 - Select the SEII PCI device driver, click remove, and confirm.

- The WinRT device category will disappear.
- Close the System Properties window.
- Reboot the computer.
- **Remove WinRT.** To do this you must know the location of the **systemroot** folder. This is on the boot drive and should be called "C:\Windows," but may be called something else depending on your installation.
 - Open **systemroot**\system\
 - Delete the **wrtdev0.vxd** file.
 - Open the **systemroot**\inf\Other folder and locate the 4pi .inf file. It should have the string "4pi" in its filename and likely will end with "FPXDRVR.INF." When this .inf file pertaining to 4pi is found, delete it.
 - Use the Run command in the START menu to run **regedit**
 - Navigate the registry tree to
HKEY_LOCAL_MACHINE→System→CurrentControlSet→Services→Class→WinRT
 - Delete the **WinRT** folder.
 - Close **regedit**.
- Proceed to the [4pi FPXDriver installation procedure](#) appropriate for your operating system.